



METAL ACTIVITY TABLE

X7000 & X8000 SERIES TABLES ASSEMBLY INSTRUCTIONS

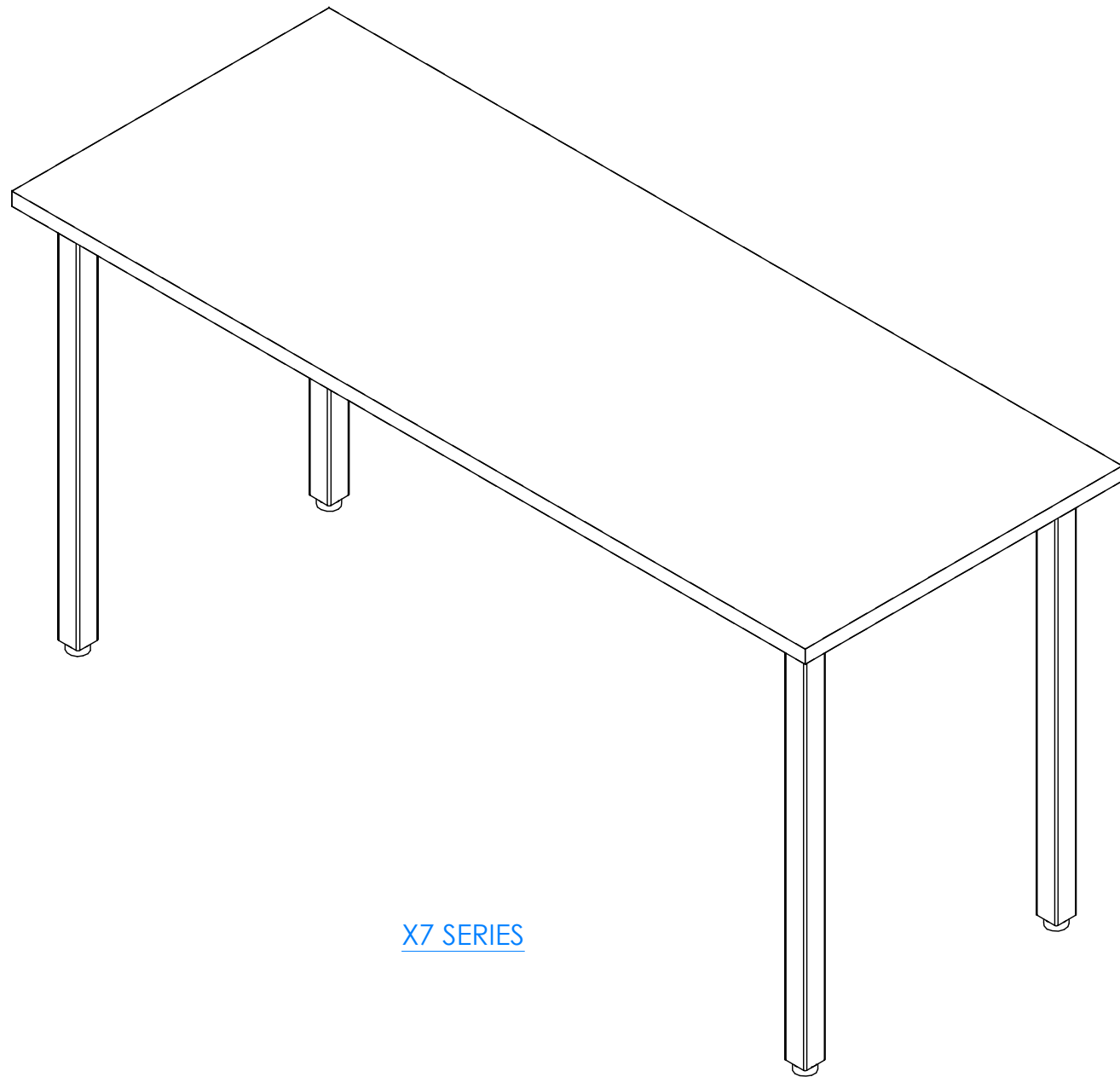

X7 SERIES

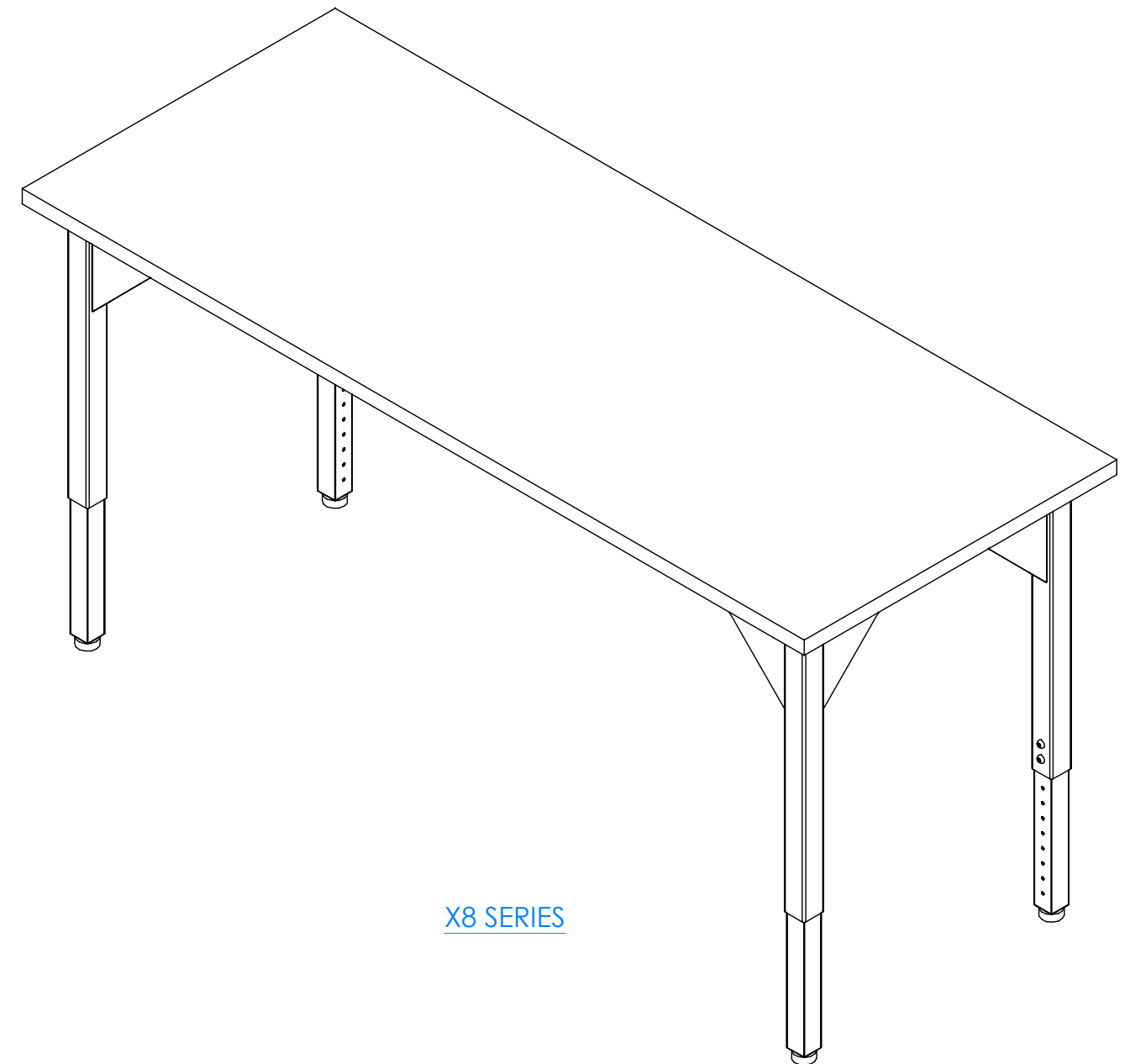

X8 SERIES

NOTE:

1. MODEL VIEWS MAY NOT REPRESENT EXACT MODEL PURCHASED

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY: EMR APR BY: APR BY:	DATE DWN: 3/3/14 DATE APR: DATE APR:
---	-----------------------------------	--

CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED.
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



TITLE: METAL ACTIVITY TABLE ASSEMBLY INSTRUCTIONS			
DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 1 OF 10	REVISION: REV B

TOOLS REQUIRED

½" WRENCH

½" SOCKET AND TORQUE WRENCH - OPTIONAL

⅛" ALLEN WRENCH

ASSEMBLY COMPONENTS

ITEMS INCLUDED	PART #	PART DESCRIPTION	QTY
TOP ASSEMBLY	VARIES	VARIES WITH STYLE	1
STEEL FRAME	VARIES	VARIES WITH STYLE	1
GUSSET - SEE NOTE	254681	GUSSET, LOBO TABLE, 4.5X4.5, SET OF 4	1
SILICONE	100711	SILICONE, SEALANT, CLEAR	1
LEG KIT - X7000 SERIES	249375	LEG KIT - FIXED HEIGHT LEGS	1
OR			
LEG KIT - X8000 SERIES	249374	LEG KIT - ADJ. HEIGHT LEGS	1

NOTE:

1. COMPONENTS VARY WITH STYLE OF TABLE PURCHASED.
 - 1.1. FRAME COLORS:
 - 1.1.1. BLACK
 - 1.2. TOP SELECTIONS:
 - 1.2.1. CHEMGUARD, PHENOLIC, 1¾" BUTCHER BLOCK MAPLE, EPOXY, SHOPTOP, OR HPL.
 - 1.3. HEIGHTS OF TABLE:
 - 1.3.1. X7 = 30"H
 - 1.3.2. X8 = ADJUSTABLE
2. GUSSETS ARE REQUIRED ON TABLES WITH PHENOLIC, EPOXY, OR BUTCHER BLOCK TOPS.
3. **XX** INDICATES ITEMS USED ONLY FOR ADJUSTABLE HEIGHT TABLES (X8000 SERIES).

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	3/3/14
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	

CONFIDENTIAL:
 ALL DESIGN RIGHTS RESERVED.
 THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



DIVERSIFIED SPACES™

TITLE:
METAL ACTIVITY TABLE TOOLS AND ASSEMBLY COMPONENTS


DRAWING NUMBER:	SCALE:	SHEET:	REVISION:
X7000 & X8000 SERIES TABLES	NTS	2 OF 10	REV B

STEP 1:

- a. IF YOU DO NOT HAVE A PHENOLIC OR EPOXY TOP, SKIP TO STEP 3.
- b. APPLY SILICONE ADHESIVE TO TOP EDGES OF METAL FRAME.
- c. APPLY SILICONE ADHESIVE IN DOT PATTERN LIKE SHOWN BELOW FOR BEST RESULTS.
- d. PLACE TOP ON METAL FRAME WITH PROPER OVERHANG (CENTERED ON FRAME).
- e. CLAMP ON TOP TO METAL FRAME TO REMOVE ANY WARP USING SHIMS AND C-CLAMPS, IF NEEDED, WHEN ALIGNING THE TOP (DO THIS BEFORE ADHESIVE SETS UP).
- f. LET CURE FOR 6-12 HOURS.
- g. **DO NOT USE SCREWS WITH PHENOLIC OR EPOXY TOPS!**

HARDWARE REQUIRED FOR STEP

QTY	DESCRIPTION	PART NUMBER
1	STEEL FRAME	VARIES
1	SILICONE	100711

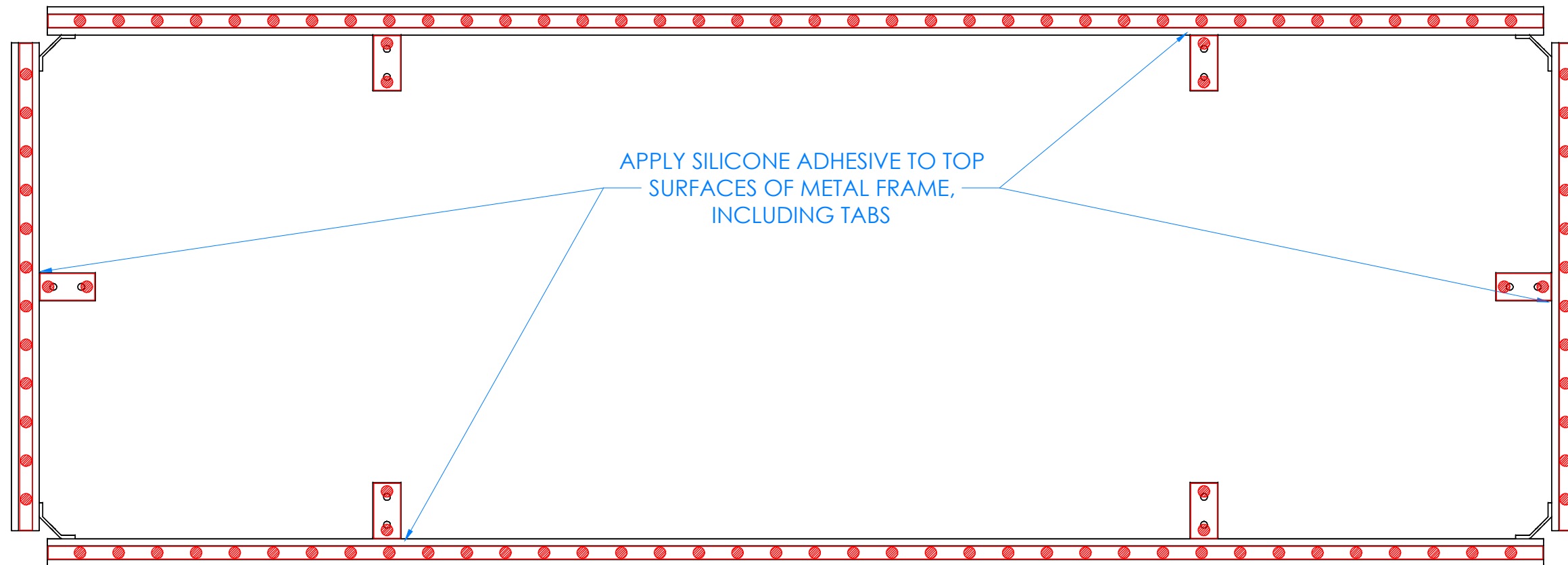

TOP OF TABLE FRAME SHOWN

NOTE:

1.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	3/3/14
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	




CONFIDENTIAL:
 ALL DESIGN RIGHTS RESERVED.
 THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

TITLE:
METAL ACTIVITY TABLE ASSEMBLY INSTRUCTIONS

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.

DRAWING NUMBER:	SCALE:	SHEET:	REVISION:
X7000 & X8000 SERIES TABLES	NTS	3 OF 10	REV B

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

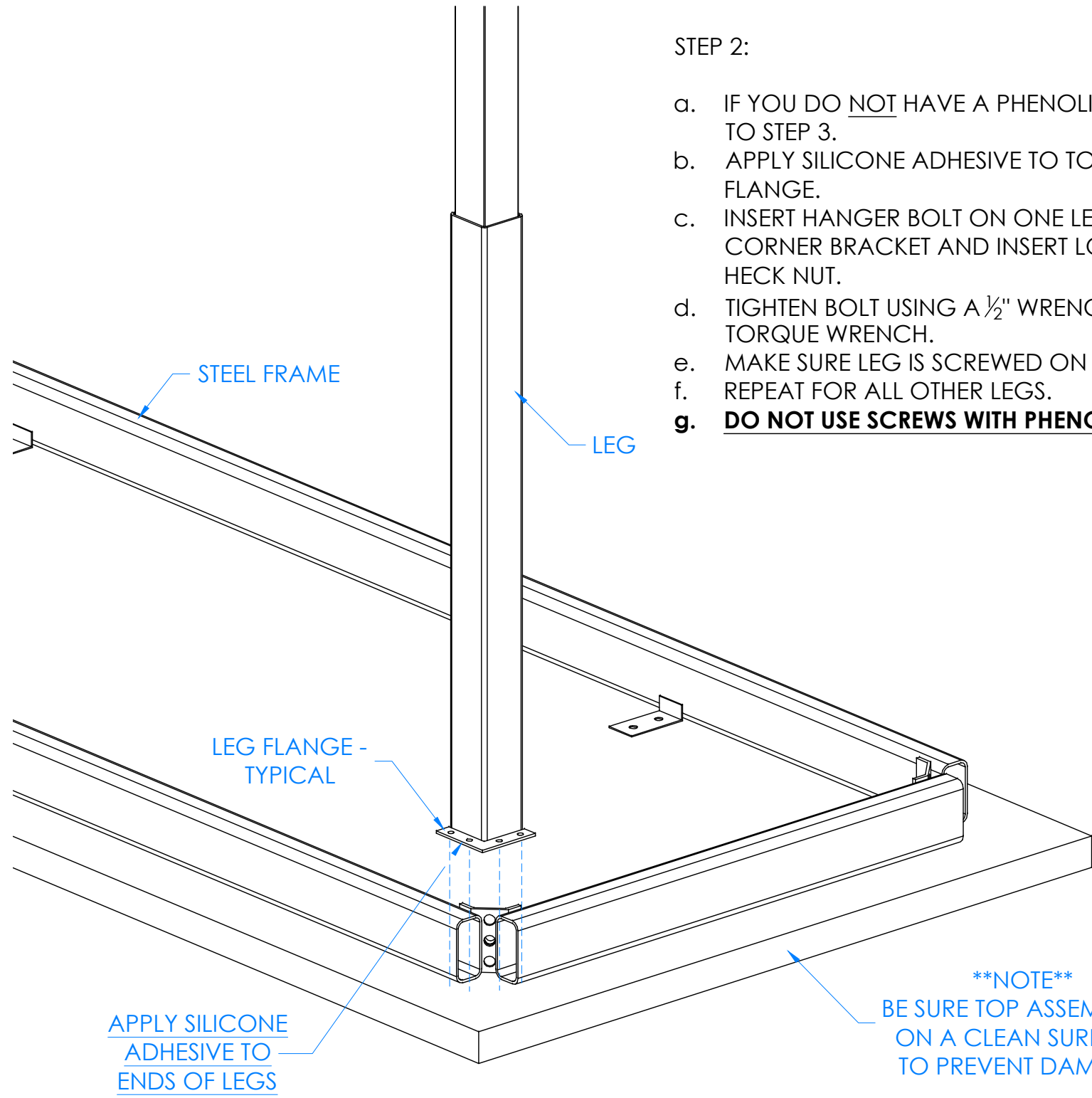


STEP 2:

- a. IF YOU DO NOT HAVE A PHENOLIC OR EPOXY TOP, SKIP TO STEP 3.
- b. APPLY SILICONE ADHESIVE TO TOP EDGES OF LEG FLANGE.
- c. INSERT HANGER BOLT ON ONE LEG THROUGH THE CORNER BRACKET AND INSERT LOCK WASHER AND HECK NUT.
- d. TIGHTEN BOLT USING A 1/2" WRENCH OR 1/2" SOCKET AND TORQUE WRENCH.
- e. MAKE SURE LEG IS SCREWED ON TIGHT!
- f. REPEAT FOR ALL OTHER LEGS.
- g. **DO NOT USE SCREWS WITH PHENOLIC OR EPOXY TOPS!**

HARDWARE REQUIRED FOR STEP

QTY	DESCRIPTION	PART NUMBER
1	LEG KIT - X7000 SERIES	249375
1	LEG KIT - X8000 SERIES	249374
1	SILICONE	100711

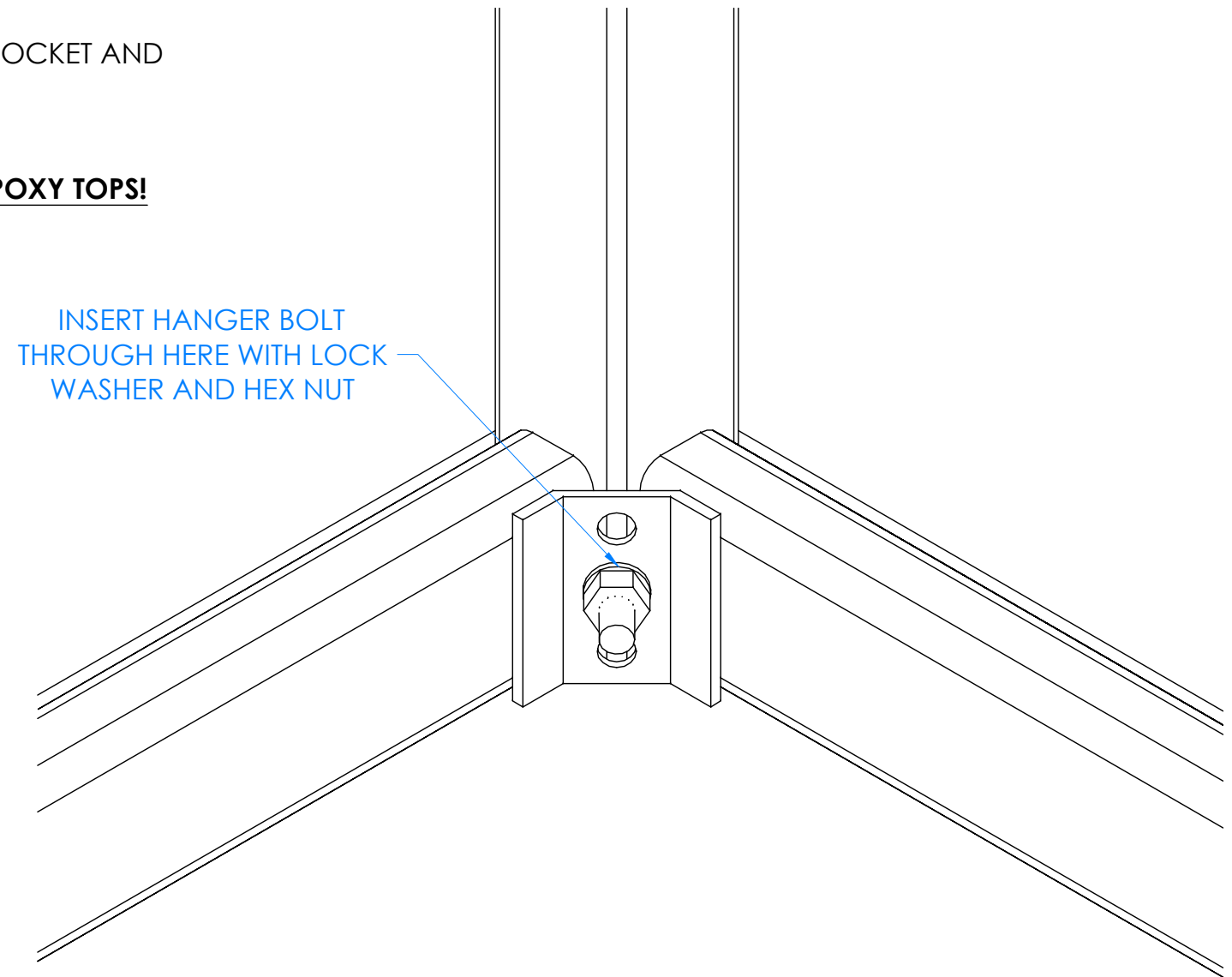

INSERT HANGER BOLT THROUGH HERE WITH LOCK WASHER AND HEX NUT

CLOSE UP OF CORNER LEG ASSEMBLY

NOTE:

1.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY: EMR APR BY: APR BY:	DATE DWN: 3/3/14 DATE APR: DATE APR:
---	-----------------------------------	--



CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED.
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

TITLE:
METAL ACTIVITY TABLE
ASSEMBLY INSTRUCTIONS

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.

DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 4 OF 10	REVISION: REV B
---	---------------	-------------------	--------------------


REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

HARDWARE REQUIRED FOR STEP

QTY	DESCRIPTION	PART NUMBER
1	TOP ASSEMBLY	VARIES

STEP 3:

- a. IF YOU HAVE A PHENOLIC OR EPOXY TOP, REFER BACK TO STEP 1 & 2
- **DO NOT USE SCREWS WITH PHENOLIC OR EPOXY TOPS!**
- b. WITH PROPER OVERHANG (CENTERED ON TOP), ATTACH FRAME TO TOP.
- c. IF YOU HAVE AN HPL, CHEMGUARD, MAPLE, OR SHOPTOP, PRE-DRILL HOLES $\frac{7}{8}$ " DEEP IN TOP USING $\frac{1}{8}$ " DRILL BIT TO PREVENT CRACKING.
- d. PLEASE MEASURE PROPER DEPTH.

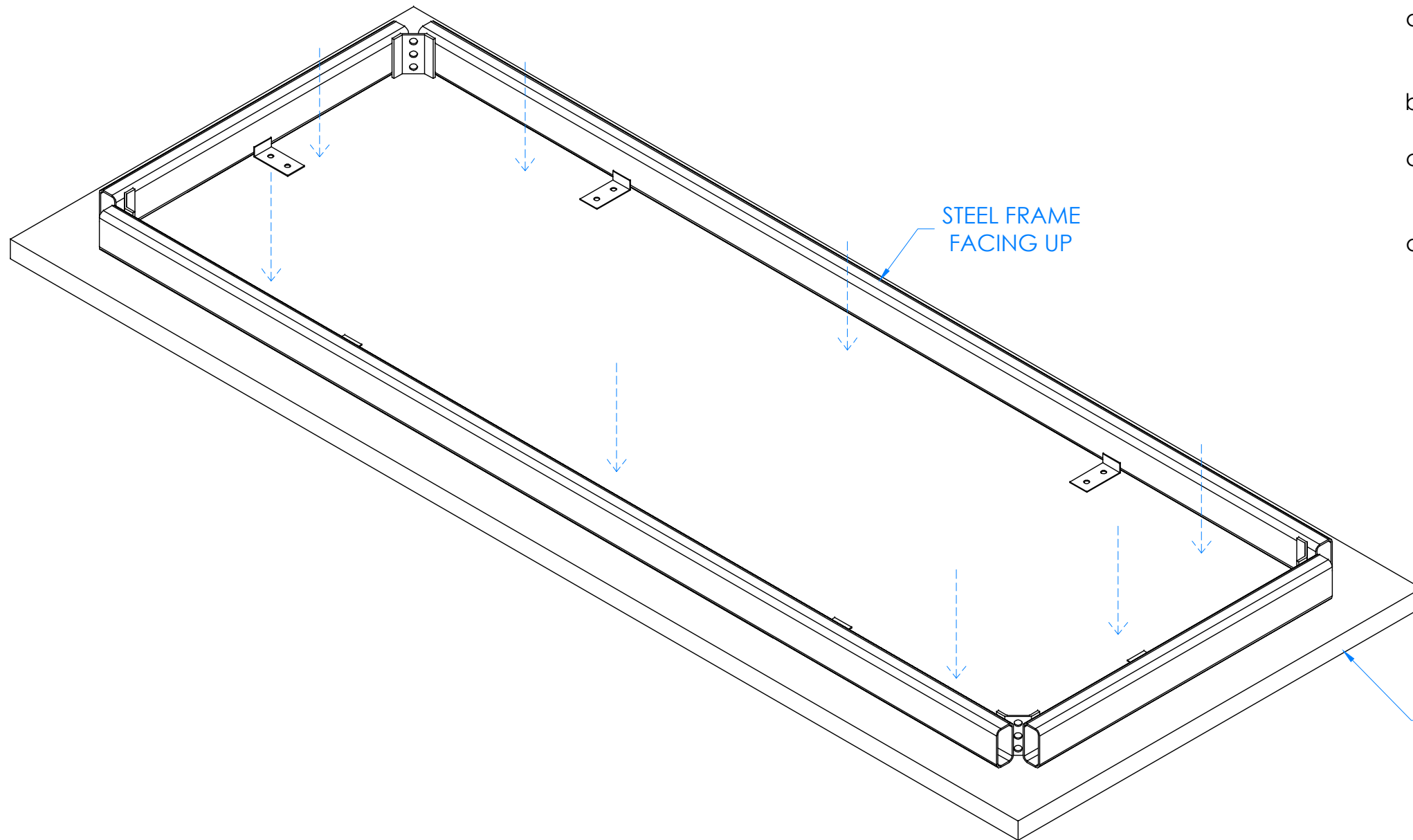

NOTE:

1.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY: EMR APR BY: APR BY:	DATE DWN: 3/3/14 DATE APR: DATE APR:
---	-----------------------------------	--



CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED.
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

TITLE:
METAL ACTIVITY TABLE
ASSEMBLY INSTRUCTIONS

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.

DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 5 OF 10	REVISION: REV B
--	---------------	-------------------	--------------------

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				


STEP 4:

- a. IF YOU HAVE A PHENOLIC OR EPOXY TOP, REFER BACK TO STEP 1 & 2
 - **DO NOT USE SCREWS WITH PHENOLIC OR EPOXY TOPS!**
- b. INSERT HANGER BOLT ON ONE LEG THROUGH THE CORNER BRACKET AND INSERT LOCK WASHER AND HEX NUT.
- c. TIGHTEN BOLT USING 1/2" WRENCH OR 1/2" SOCKET AND TORQUE WRENCH.
- d. SCREW LEGS TO TOP BY ATTACHING WITH HARDWARE PROVIDED AND SCREWING THRU LEG FLANGE INTO TOP.
- e. PRE-DRILL HOLES TO PREVENT CRACKING.
- f. MAKE SURE LEG IS SCREWED ON TIGHT.
- g. REPEAT FOR OTHER LEGS.


HARDWARE REQUIRED FOR STEP

QTY	DESCRIPTION	PART NUMBER
1	LEG KIT - X7000 SERIES	249375
1	LEG KIT - X8000 SERIES	249374

INSERT HANGER BOLT THROUGH HERE WITH LOCK WASHER AND HEX NUT

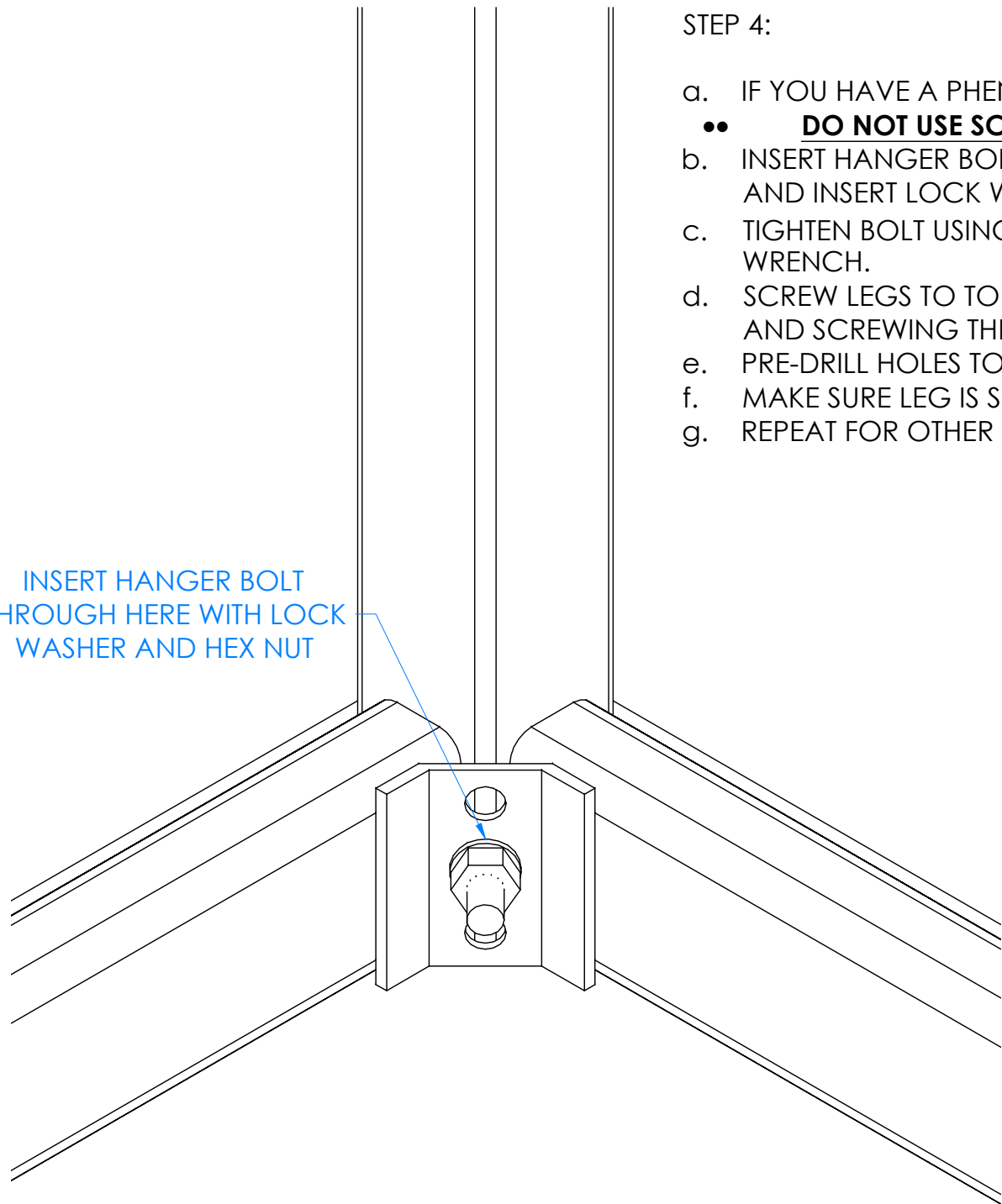

CLOSE UP OF CORNER LEG ASSEMBLY

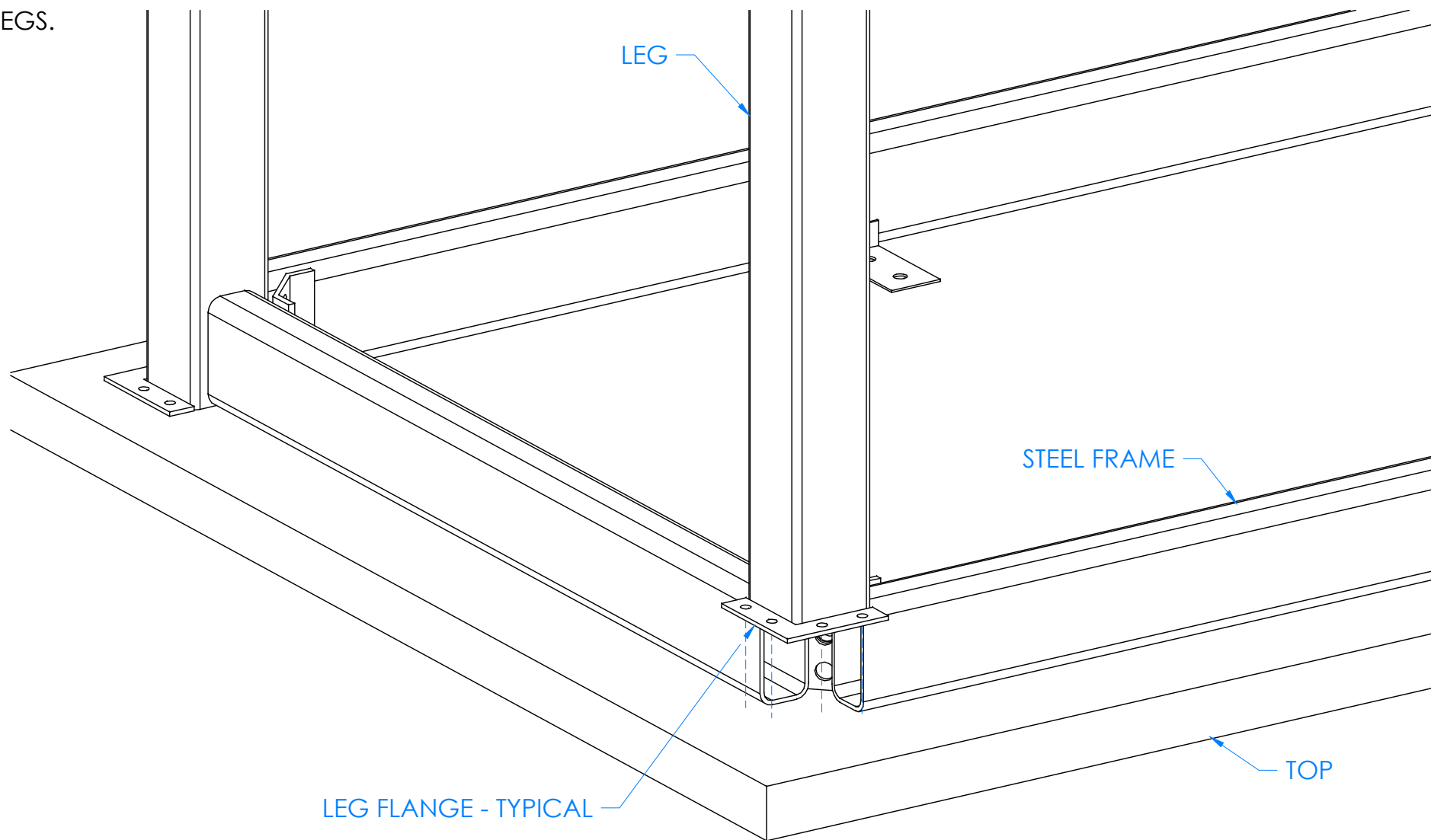

LEG FLANGE - TYPICAL

NOTE:

1.

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE	DWN BY: EMR	DATE DWN: 3/3/14
DECIMALS: .X = ± .030	APR BY:	DATE APR:
.XX = ± .020	APR BY:	DATE APR:
.XXX = ± .010	APR BY:	DATE APR:
ANGLES: = ± 1°		
CABINET SQUARENESS = ± .0625		




CONFIDENTIAL: ALL DESIGN RIGHTS RESERVED. THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

TITLE: METAL ACTIVITY TABLE ASSEMBLY INSTRUCTIONS

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.

DRAWING NUMBER:	SCALE:	SHEET:	REVISION:
X7000 & X8000 SERIES TABLES	NTS	6 OF 10	REV B

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

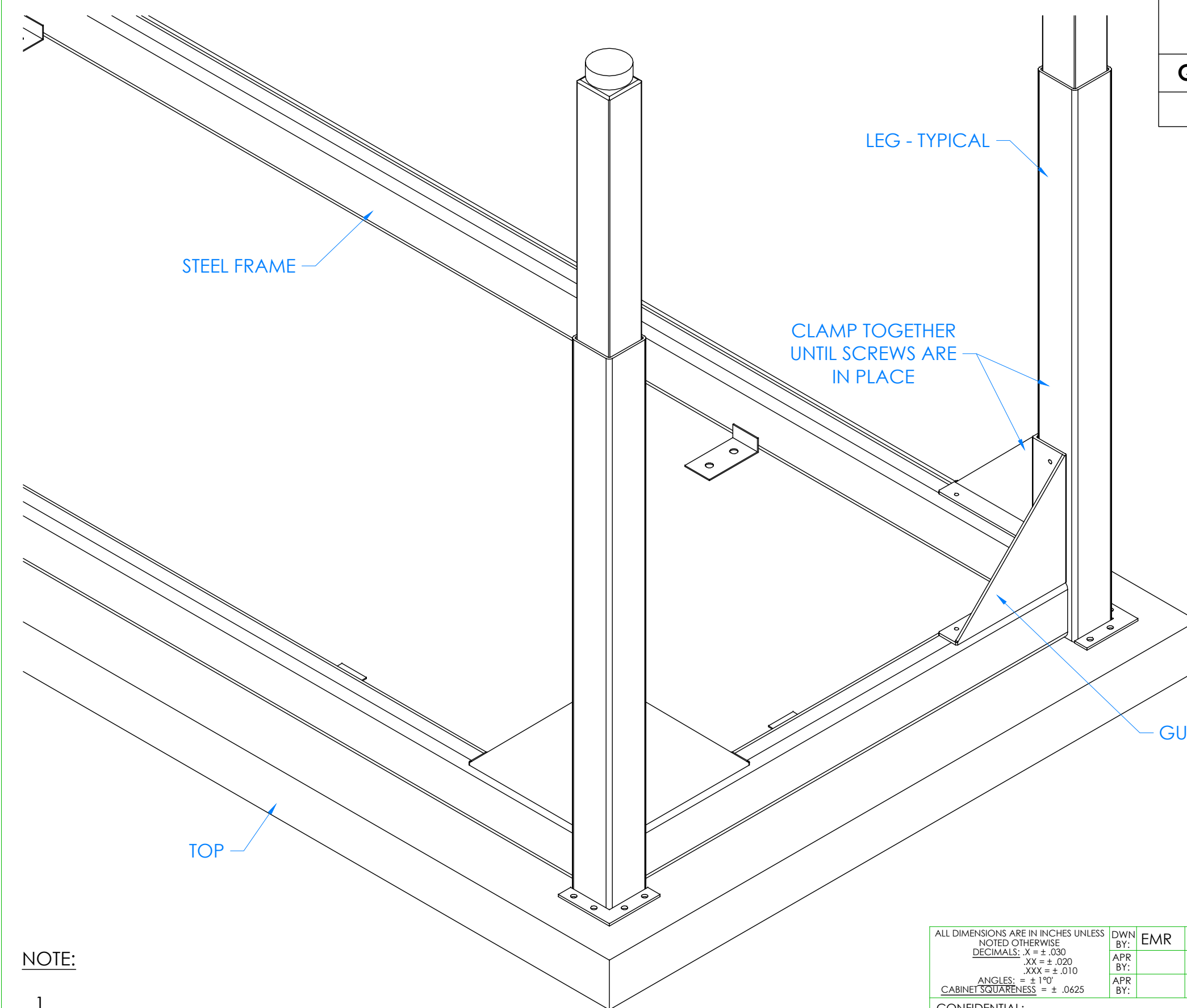

HARDWARE REQUIRED FOR STEP

QTY	DESCRIPTION	PART NUMBER
1	GUSSET - SET OF 4	254681

STEP 5:

- a. IF YOUR TABLE WAS SUPPLIED WITH LEG GUSSETS, PLACE A GUSSET IN EACH CORNER AROUND EACH LEG.
- b. CLAMP ONE SIDE OF GUSSET TO LEG.
- c. ATTACH GUSSET INTO PRE-DRILLED HOLES ON STEEL FRAME WITH HARDWARE PROVIDED.
- d. REPEAT FOR REMAINING GUSSETS.

NOTE:
1.

CLOSE UP OF GUSSET ASSEMBLY SHOWN

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				


ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY: EMR APR BY: APR BY:	DATE DWN: 3/3/14 DATE APR: DATE APR:
---	-----------------------------------	--

CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.
THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



TITLE:
METAL ACTIVITY TABLE
ASSEMBLY INSTRUCTIONS

DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 7 OF 10	REVISION: REV B
--	---------------	-------------------	--------------------

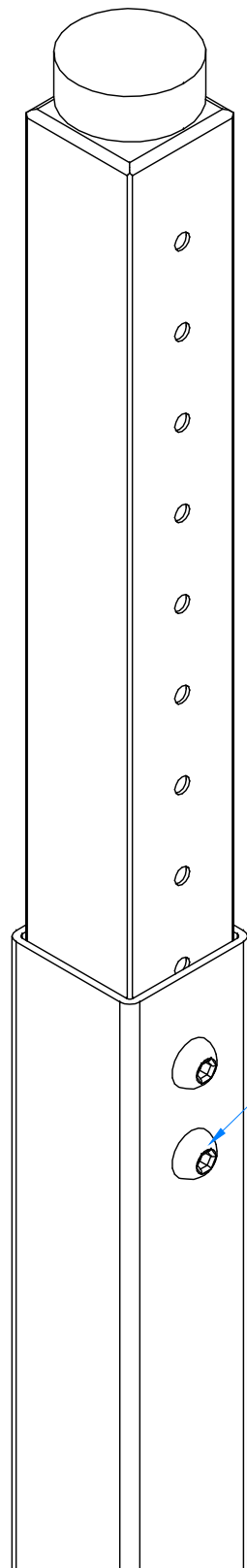

STEP 6:

- a. IF YOU DO NOT HAVE THE X8000 SERIES, SKIP TO STEP 7.
- b. ADJUST LEGS TO DESIRED HEIGHT BY REMOVING TWO ALLEN SCREWS IN EACH LEG.
- c. BE CAREFUL TO SUPPORT THE LEG AS IT WILL FALL ONCE SCREWS ARE REMOVED.
- d. MOVE LEG TO A NEW HEIGHT AND RE-INSERT ALLEN SCREWS AND TIGHTEN.
- e. REPEAT FOR OTHER LEGS.

ALLEN SCREW - TYPICAL

NOTE:

1.

CLOSE UP OF X8000 SERIES LEG ASSEMBLY

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY:	EMR	DATE DWN:	3/3/14
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	
	APR BY:		DATE APR:	


CONFIDENTIAL:
 ALL DESIGN RIGHTS RESERVED
 THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.
 THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



DIVERSIFIED SPACES™

TITLE:
METAL ACTIVITY TABLE ASSEMBLY INSTRUCTIONS

DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 8 OF 10	REVISION: REV B
---	---------------	-------------------	--------------------

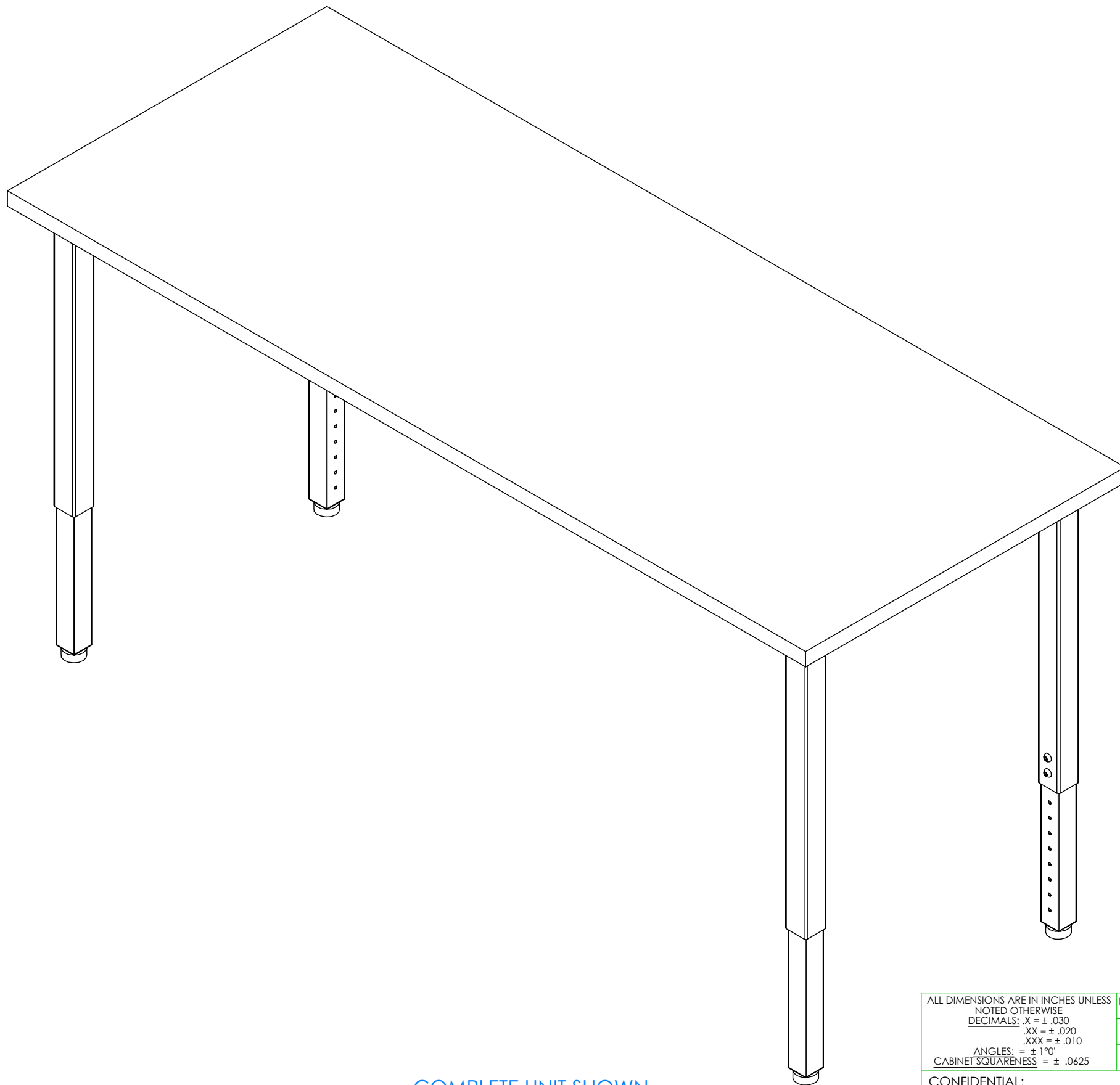

STEP 7:

- a. WITH HELP FROM ANOTHER PERSON, TURN TABLE OVER SO TOP IS UP.
- b. SET YOUR TABLE IN DESIRED LOCATION.
- c. ADJUST LEG LEVELERS TO LEVEL TABLE.

NOTE:

1.

COMPLETE UNIT SHOWN

REV.	DESCRIPTION	DATE	NAME	ECN No.
B	UPDATED SILICONE PROCEDURE	8/31/23	RJM	N/A
A	CREATED DRAWING	3/3/14	EMR	N/A
REVISION LEVEL				

ALL DIMENSIONS ARE IN INCHES UNLESS NOTED OTHERWISE DECIMALS: .X = ± .030 .XX = ± .020 .XXX = ± .010 ANGLES: = ± 1° CABINET SQUARENESS = ± .0625	DWN BY: EMR APR BY: APR BY:	DATE DWN: 3/3/14 DATE APR: DATE APR:
---	-----------------------------------	--

CONFIDENTIAL:
ALL DESIGN RIGHTS RESERVED
THE INFORMATION, DATA, DESIGN, AND DRAWINGS EMBODIED IN THIS DOCUMENT AND ALL REPRODUCTIONS THEREOF ARE THE EXCLUSIVE PROPERTY OF DIVERSIFIED WOODCRAFTS, INC.

THESE DOCUMENTS ARE SUPPLIED WITH THE UNDERSTANDING THAT THEY WILL NOT BE DISCLOSED TO THIRD PARTIES WITHOUT WRITTEN CONSENT OF DIVERSIFIED WOODCRAFTS, INC. ANY REPRODUCTION IN PART OR WHOLE WITHOUT WRITTEN CONSENT IS SUBJECT TO LEGAL ACTION.



TITLE: METAL ACTIVITY TABLE ASSEMBLY INSTRUCTIONS		DRAWING NUMBER: X7000 & X8000 SERIES TABLES	SCALE: NTS	SHEET: 9 OF 10	REVISION: REV B
---	--	--	---------------	-------------------	--------------------